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SAFETY, STORAGE AND SETUP INFORMATION

1) **General Information:** the following general precautions for handling must be observed during all phases of operation and service of CAEN equipment. Failure to comply with these precautions or with specific warnings elsewhere in the product manual could result in personal injury or damage to the equipment. The precautions listed below represent warnings of certain dangers of which CAEN is aware. The user of the product, should follow these warnings and all other safety precautions necessary for the safe operation of the equipment in the operating environment.

CAEN shall not be liable for user’s failure to comply with these requirements

2) **Conventions Used in This paragraph:**

CAUTION indicates a hazard that could damage the equipment. Such damage may invalidate the warranty.

3) **Safety Terms and Symbols on the Product:** these terms may appear on the product:

- **DANGER:** indicates a hazardous situation which, if not avoided, could result in serious injury or death.
- **WARNING:** indicates a hazardous situation which, if not avoided, could result in death or serious injury.
- **CAUTION:** indicates a situation or condition that, if not avoided, could cause physical injury or damage the product and/or its environment.

The following symbols may appear on the product:

 Caution, refers to the product manual	 Caution, risk of electric shock	 Productive Conductor Terminal
 Earth (Ground) Terminal	 Alternating Current	 Three-Phase Alternating Current

4) **Unpacking and Inspection:**

CAUTION: observe precautions for handling electrostatic sensitive devices.

- Inspect containers for damage during shipment. Report any damage to the freight carrier for possible insurance claims.
- Check that all the components received match those listed on the enclosed packing list. (CAEN cannot accept responsibility for missing items unless we are notified promptly of any discrepancies.)
- Open shipping containers; be careful not to damage any contents.
- Inspect contents and report any damage. The inspection should confirm that there is no exterior damage to the unit such as broken knobs or connectors and that the front panel and display face are not scratched or cracked. Keep all packing material until the inspection has been completed.
- If damage is detected, file a claim with carrier immediately and notify CAEN service.
- If equipment must be returned for any reason, carefully repack equipment in the original shipping container with original packing materials if possible. Please contact CAEN service.

- If equipment is to be installed later, place equipment in original shipping container and store in a safe place until ready to install.

5) **ESD precautions:** CAEN equipment may be classified as ESD sensitive item¹.

CAUTION: during long term or temporary storage the items shall be covered by ESD protective bags. Handling of the items without ESD protective covering shall be performed only into approved ESD Protected Area (EPAs); any operator shall be properly grounded (e.g., by wrist-straps) before handling the item. Avoid handling operations if the relative humidity is below 30% or take extra ESD safety precautions (such as redundant ground clips) if such handling is unavoidable.

¹ Excluded all the devices provided with independent case or chassis. Devices that consist of a circuit board with no case or chassis (e.g. VME, CAMAC, NIM modules or HV/LV Power supply boards) comply with the EMC directive if installed in a CE marked system.

6) **Storage Condition²:** CAEN equipment shall be handled and stored in a clean, dust free environment under the following conditions:

	Minimum	Maximum
Temperature	- 30° C	+ 80° C
RH (non-condensing)	5 %	90 %

²Typical ranges, for more details see the product manual.

7) **Handling:** CAEN equipment shall be handled only with lint-free gloves (e.g. nitrile); never touch the item with bare, ungloved hands.

CAUTION: do not subject the item to undue shock of vibrations; do not bump, drop or slide shipping containers. Do not leave items or shipping containers unsupervised in areas where untrained personnel may mishandle the items.

8) **Instruction for Cleaning:** if required, the equipment may be cleaned with isopropyl alcohol or deionised water and air dried. Clean the exterior of the product only. Do not apply cleaner directly to the items or allow liquids to enter or spill on the product.

Cleaning the Touchscreen (if present)

In order to clean the touchscreen, wipe the screen with a towelette designed for cleaning monitors or with a clean cloth moistened with water. Do not use sprays or aerosols directly on the screen; the liquid may seep into the housing and damage a component. Never use solvents or flammable liquids on the screen.

Cleaning the air vents (if present)

It is recommended to clean the air vents (if present) occasionally on all vented sides of the board. Lint, dust, and other foreign matter can block the vents and limit the airflow. Be sure to unplug the board before cleaning the air vents and follow the general cleaning safety precautions.

CAUTION: CAEN recommends cleaning the device using the following precautions:

- Never use solvents or flammable solutions to clean the board.
- Never immerse any parts in water or cleaning solutions; apply any liquids to a clean cloth and then use the cloth on the component.
- Always unplug the board when cleaning with liquids or damp cloths.
- Always unplug the board before cleaning the air vents.
- Wear safety glasses equipped with side shields when cleaning the board

9) **Transport:**

CAUTION: CAEN equipment can be safely transported in their original shipping packages.

10) Installation and Operation: only qualified personnel should perform installation, operation, and maintenance procedure.

11) Safety Precautions: keep extreme caution when an electrical shock hazard is present. Lethal voltages may be present on cable connector jacks or test fixtures. Warning label indicates HV sources which are hazardous to personnel; during test or integration all exposed HV connections shall be insulated or protected.

Before operating an instrument, make sure the line cord is connected to a properly grounded power receptacle. Inspect the connecting cables, test leads, and jumpers for possible wear, cracks, or breaks before each use.

When installing equipment, where access to the main power cord is restricted, such as rack mounting, a separate main input power disconnect device must be provided, in close proximity to the equipment and within easy reach of the operator. It is recommended to disconnect ac line from the entire test system and discharge any capacitors before:

- connecting or disconnecting cables or jumpers
- installing or removing switching cards
- removing cover and making internal changes, such as installing or removing jumpers, if required
- **CAUTION:** The instrument and accessories must be used in accordance with its specifications and operating instructions, or the safety of the equipment may be impaired.
- **CAUTION:** Do not exceed the maximum signal levels of the instruments and accessories, as defined in the specifications and operating information.
- **CAUTION:** When fuses are used in a product, replace with same type and rating for continued protection against fire hazard.

12) Ventilation Requirements

CAUTION: all power supplied CAEN NIM, VME and CAMAC modules must be used only with crates with forced cooling air flow.

CAEN modules without power supply, such as Attenuators and Delay units, do not require crates with forced cooling air flow.

CAUTION: rack bound CAEN HV Power Supply Systems must be installed with a minimum 10 cm space both above and below the system.

13) Medical Use: CAEN equipment is not designed with components and testing procedures that would ensure a level of reliability suitable for use in treatment and diagnosis of humans. These items are not intended to be a substitute for any form of established process or equipment used to monitor or safeguard human health and safety in medical treatment.

14) Operating Environmental Requirements²

	Minimum	Maximum
Temperature	0° C	+ 40° C
RH (non-condensing)	10 %	90 %
Environmental	Indoor use	
Altitude	<2000m	
Pollution degree	2	
Overvoltage category	II	
EMC environment	Commercial and light industrial	
IP degree	IPX0, not for wet location	

² Typical ranges, for more details see the product manual.

15) Installing Devices with No Case or Chassis: (e.g., VME, CAMAC, NIM modules or HV/LV Power supply boards):

CAUTION: CAEN strongly recommends to use an antistatic wrist strap and a conductive foam pad when installing these items. After removing the devices from the system or its protective antistatic wrapper, place the component flat on a grounded, static-free surface (and in the case of a board, component side up). Do not slide the component over any surface. Do not stack units on top of each other.

CAUTION: crate power should be turned off during insertion or removal of modules³ to avoid possible damage caused by **momentary** misalignment of contacts.

CAUTION: the safety of any system that incorporates the device falls under the responsibility of the assembler of the system

³ Excluded modules supporting live insertion/extraction into/from the crate.

CAUTION: to insert a module in the relative crate follow these steps:

- Choose a convenient I/O station.
- Hold the board firmly with both hands (if HV/LV Power supply boards).
- Align the PCB back side with one crate free slot rails.
- Start sliding gently the board along the rails.
- Pay attention to closer boards: insert the module carefully to avoid contacting boards to the right or left.
- Ensure that correct alignment is kept.
- Continue sliding the board holding it firmly.
- Ensure that sliding is smooth.
- Plug the board into Crate backplane connectors with gentle finger pressure.
- Ensure the correct backplane connectors mating.
- Complete installation by screw-locking the board front panel on the crate.

16) Repairs: may be executed by CAEN service only.

17) EMI considerations: CAEN equipment has been designed in compliance with the harmonized standards concerning Electromagnetic compatibility. Devices meet requirements concerning EMC. Specific test was performed in normal use condition in order determine the effective EM effects. Adhering the following rules will help EMI

- Power supply cable routing. The power supply cable is usually susceptible to conduct EMI inside the device and coupling into power supply cord. It is recommended to keep the power supply cord far from SIP/SOP cables and fixed in place with cable ties, if necessary
- SIP/SOP connector shield (if available). Sometimes the I/O connector realize poor contact between metal housing and connector plate. It is recommended to check electrical continuity
- Minimize the differential mode loop of signal wires by keeping the wires close together. This will minimize the magnetic field and reduce radiated EMI.
- Minimize the common mode capacitance of wires by keeping the length as short as possible
- If specified, connect the chassis earth connection properly
- If applicable, the devices are designed to run with spread spectrum clocking enabled: it is recommended because often produces low EM emissions.
- If not necessary, disable unused clocks. Clock signals that have no load can have high level of ringing that can lead to EMI problems. For the same reason it is recommended to disable unused HV channels or input acquisition channels

CAUTION: If the customers modify the hardware/software/firmware factory configuration, the EMC compliance will not be guaranteed.

18) Device Decommissioning: After its intended service, it is recommended to perform the following actions:

- Detach all the signal/input/output cable
- Wrap the device in its protective packaging
- Insert the device in its packaging (if present)

CAUTION: The device shall be stored only at the declared environment condition, otherwise performance and safety will not be guaranteed.

19) Disposal: The disposal of the equipment must be managed in accordance with Directive 2012/19 / EU on waste electrical and electronic equipment (WEEE).

The crossed bin symbol indicates that the device shall not be disposed with regular residual waste.

PRODUCT SUPPORT SERVICE AND REPAIR

For technical assistance on hardware and software, please follow the following procedure


- 1) **MyCAEN Account:** A MyCAEN account is required for ask support about CAEN products and register a firmware/software license. Making an account is totally free and it is possible to upgrade to MyCAEN+ anytime.

New user registration

- Go to registration page and fill the registration form with your data.
- Validate your email address by clicking on the confirm link and set a password.
- Log in to CAEN website to activate your account.

Do I have a MyCAEN+ account?

If you are already registered and need to know if you have a MyCAEN+ accounts, you will be able to check it by logging in to CAEN website, go to your dashboard and see the title of the section that could be MyCAEN or MyCAEN+.

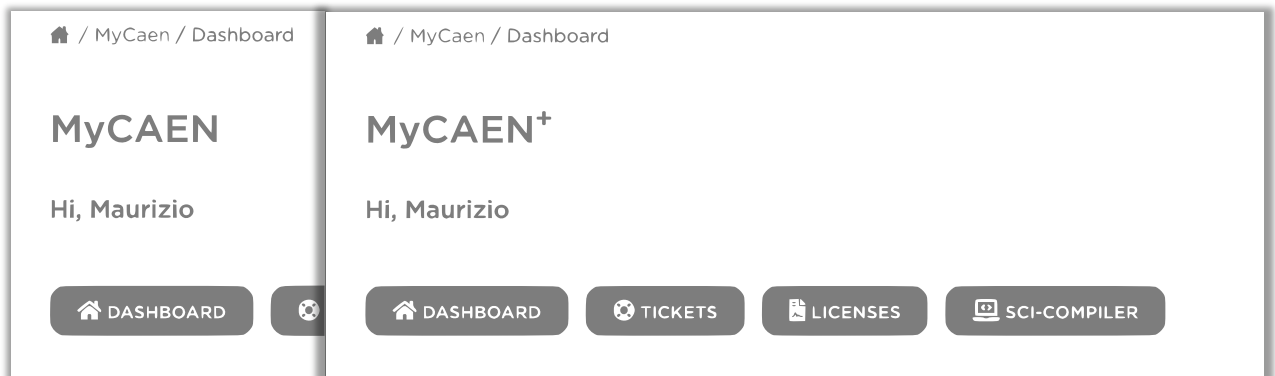Upgrade to MyCAEN+

If you are already registered on CAEN website and you have a MyCAEN account, you will just need to update to MyCAEN+ by following the instructions below:

- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button at the right top.

- Go to "Edit Profile" and on "Profile Information" update your Institutional email, country/region and Company/institute fields.
- Click on "Save Profile".

Known institutes are automatically certified by CAEN website. New institutes or generic email addresses need to be validated by CAEN staff. This operation can take up to 48 hours.


2) **Register a Device.** The product registration is required for access to support services.

Add a device

- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button on the top right.
- On the menu area click on “Dashboard”, then “Add Device” button


- Select an identification method, PID or S/N. Please refer to chapter “Find your S/N or PID” to discover how to find your board identification number.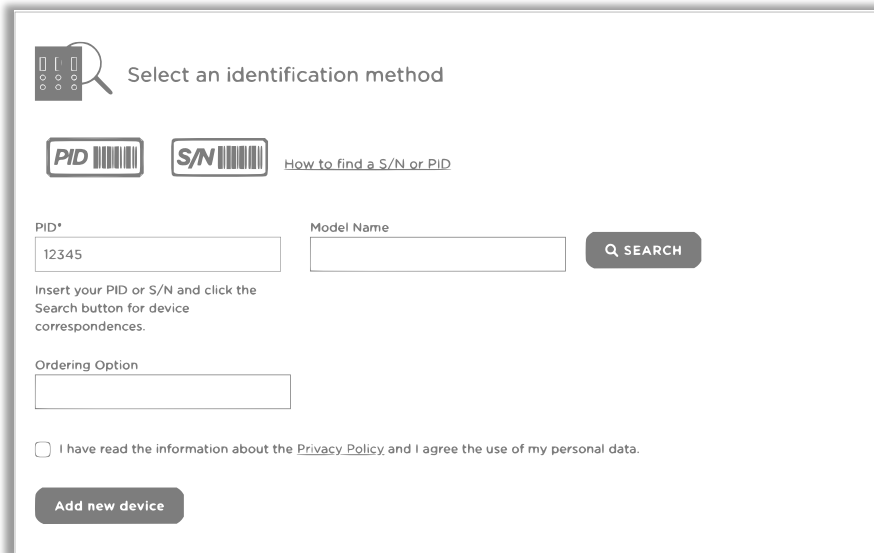


The image shows a web form titled 'Select an identification method'. It features two options: 'PID' with a barcode icon and 'S/N' with a barcode icon and a link 'How to find a S/N or PID'. Below these are input fields for 'PID*' (containing '12345') and 'Model Name'. A 'SEARCH' button is to the right. Below the input fields, there is a note: 'Insert your PID or S/N and click the Search button for device correspondences.' and an 'Ordering Option' field. At the bottom, there is a checkbox for 'I have read the information about the Privacy Policy and I agree the use of my personal data.' and an 'Add new device' button.

- Click on “Search” button and then “This is my device.”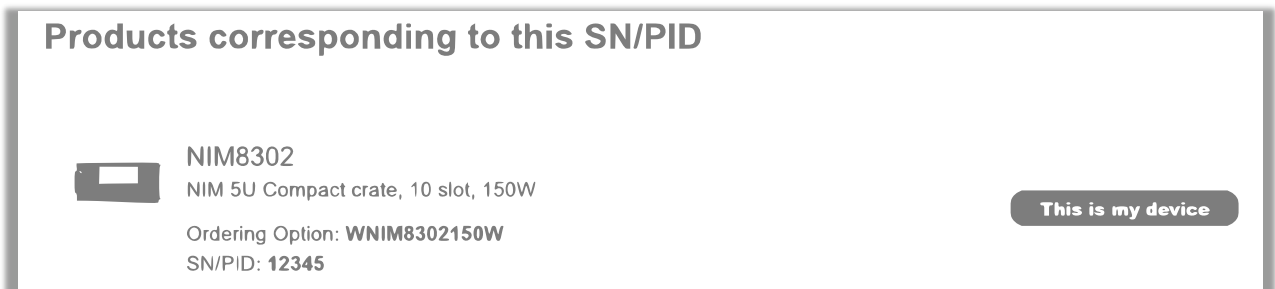

The image shows a search result titled 'Products corresponding to this SN/PID'. It displays a product image of a NIM8302 NIM 5U Compact crate, 10 slot, 150W. Below the image, the text reads: 'Ordering Option: WNIM8302150W' and 'SN/PID: 12345'. A 'This is my device' button is located to the right of the product information.

- The fields will be automatically filled by the system. If all information are correct and the Privacy Policy is accepted, it will be possible to click on “Add new device”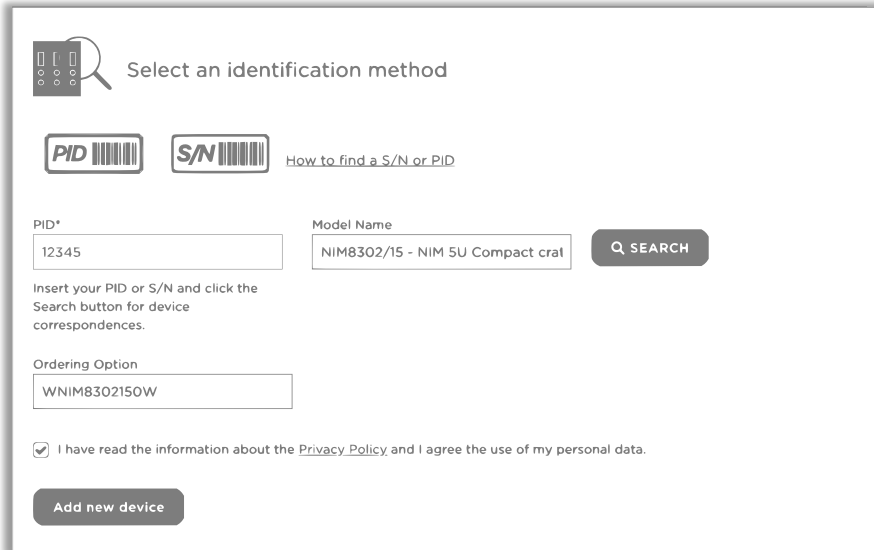

Select an identification method

PID **S/N** [How to find a S/N or PID](#)

PID* Model Name

Insert your PID or S/N and click the Search button for device correspondences.

Ordering Option


I have read the information about the [Privacy Policy](#) and I agree the use of my personal data.

If the device cannot be found, please double check your identification number; if the search gives no results instead, it will still be possible to create a support ticket without registering the device: the model name and the Serial Number or PID can be specified in the message.

Find your Serial Number (S/N) or PID

The product information label contains your Serial Number (S/N) or Product IDentifier (PID). Serial Number can be composed by 2 or more numbers, PID are composed by at least 5 char. It is located either on the back of your CAEN module, in the front or under the right side of the board.


A detail view by form factor is the following.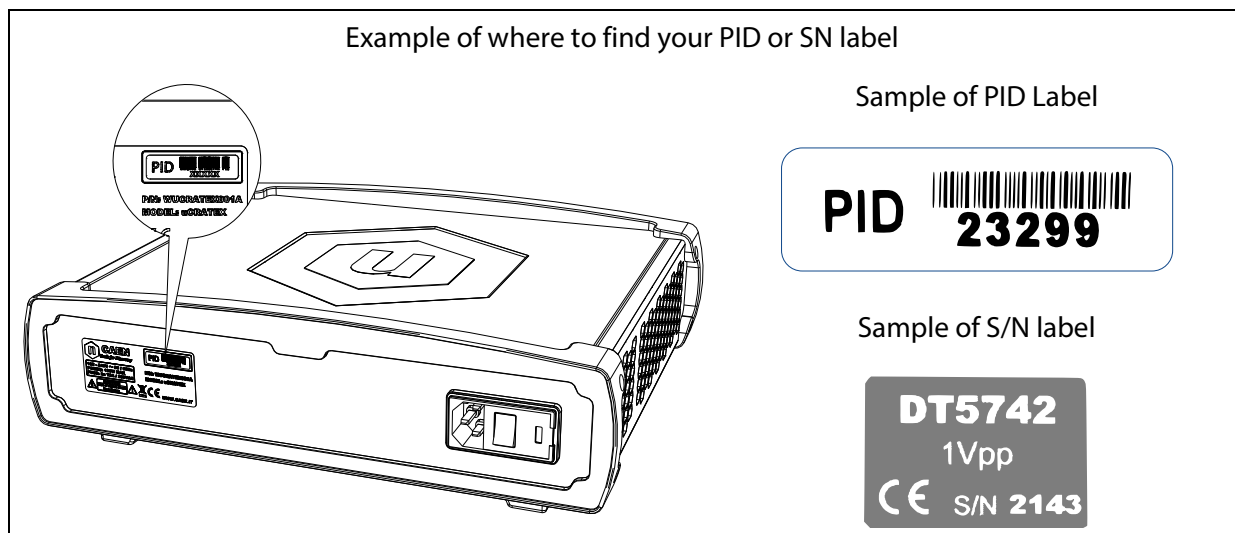

Download device list (.csv)

It is possible to download the device list in .csv format, by the following:


- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button at the right top.
- On the menu area click on "Dashboard", then "Download devices list" written on the right.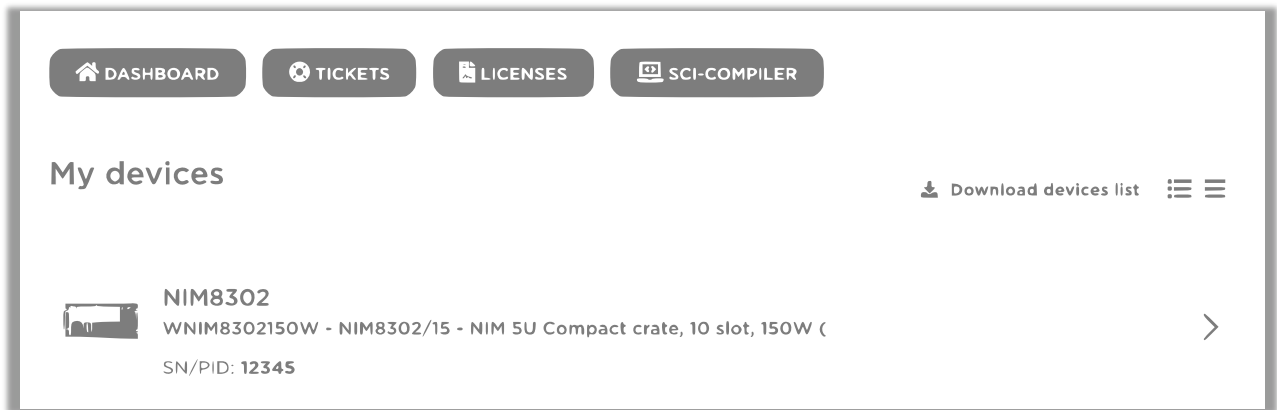

3) Support Ticketing

Create a Ticket

- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button at the right top.
- On the menu area click on "Tickets", then "Create Ticket" button.

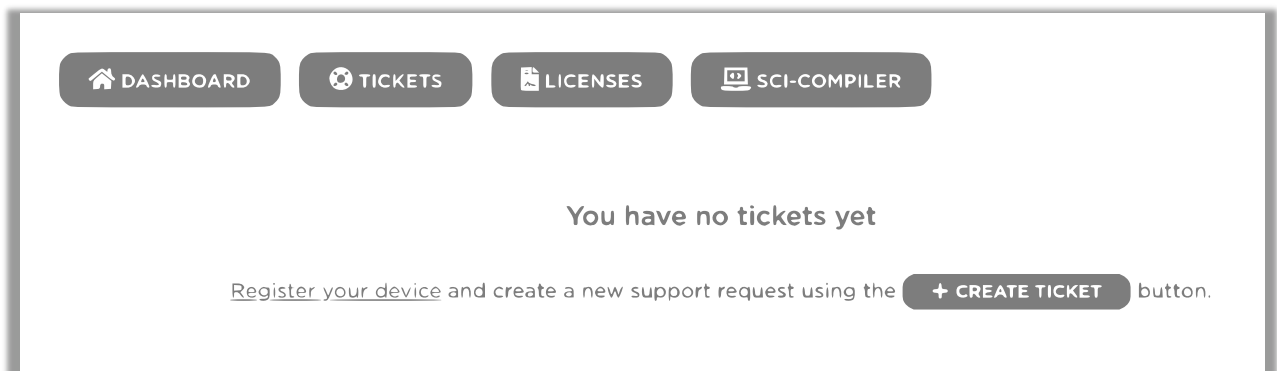- Fill all the required field being careful to select your hardware device.

When the CAEN team reply to your tickets you will receive a notification email with a direct link to the relative ticket.

Reply to a Ticket

- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button at the right top.
- On the menu area click on "Tickets", then click on the tickets you want to reply.
- Fill the form at the bottom of the page.

When the CAEN team reply to your tickets you will receive a notification email with a direct link to the relative ticket.

RMA – Return Merchandise authorization

- Create a Ticket describing your hardware failure.
- If your problem is evaluated by CAEN team as not solvable by the user, the RMA form will be activated and a notification email will be sent to the user.
- Click on the link received by email and then click on "Return Device" to fill the form with all return information.

Depending on where you are located, the product must be shipped directly to CAEN S.p.A. headquarter or to your Country Dealer. To find the proper address please visit the page at

<https://www.caen.it/sales-network/>

4) License. The user can register the software license

SCI-Compiler License

- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button at the right top.
- On the menu area click on "License", then "Activate SCI-Compiler License".
- Copy the OTP you've received by email or by mail with the USB dongle.
- If the license is valid, a new button "SCI-Compiler" will appear in the menu area.

Web License

- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button at the right top.
- On the menu area click on "License", then "Register License".
- Complete the wizard procedure by selecting the module in your possess.

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